



Resource CD Installation Help

[I. Installing USB drivers](#)


[II. Installing PolyTemp™](#)

[III. Installing TDMS Viewer](#)

I. Installing USB drivers

IMPORTANT: TO CONNECT TO THE POLYSCIENCE CIRCULATING BATH USING USB, INSTALL THE USB DRIVER BEFORE PLUGGING THE UNIT INTO YOUR PC.



- 1) Insert the PolyScience Driver CD into your PC.
- 2) Click the "Install USB Driver Icon"
- 3) A command-line window will briefly appear while the driver is installing. Installation is complete when this screen closes.

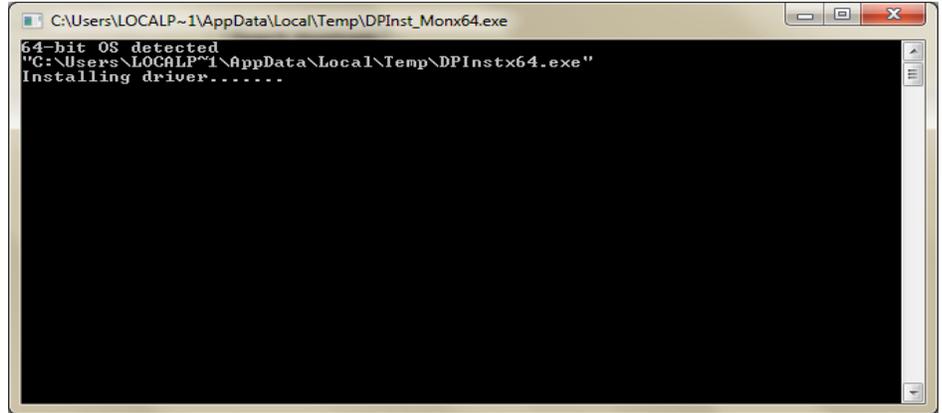

- 4) Plug the USB cable connected to the Circulating Bath into your PC. Windows will complete the driver installation.


Installation notes:

- Your circulating bath will be installed as a Virtual COM port.
- The default designation is COM3.
- If you have more than 3 physical serial ports or additional virtual COM devices, the next available COM port will be designated.
- To identify which port is assigned to the Circulating Bath, right-click on My Computer and select Properties. Click Device Manager.
- The Circulating Bath USB connection will be listed under "Ports (COM & LPT)". If you have multiple listings for COM ports, double-click on each one and check the Manufacturer on the general tab. The Circulating Bath will list "FTDI" in the manufacturer field, referring to the USB chipset type in your PolyScience unit.

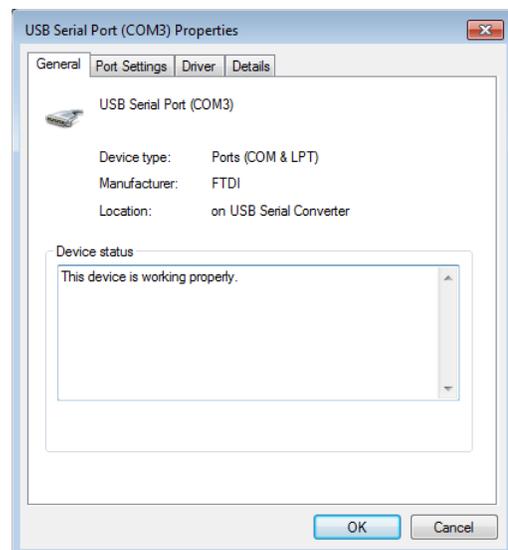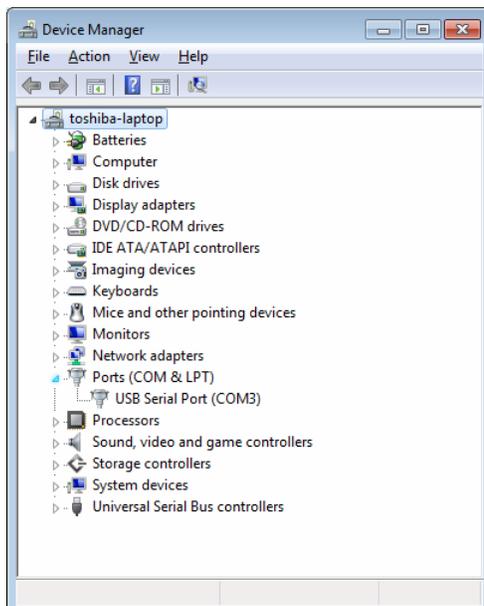


II. Installing PolyTemp™



- 1) Insert the PolyScience Driver CD into your PC.
- 2) Click the "Install PolyTemp™ Icon"
- 1) Click "Next"

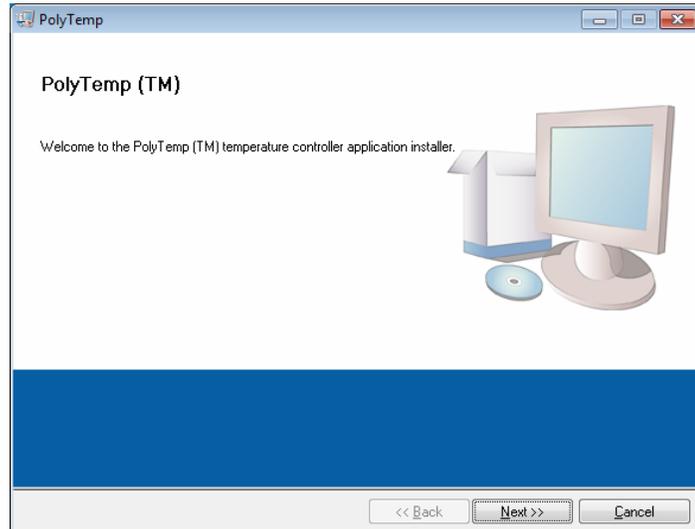


- 2) Choose the directories that the installation files will be installed in. It is recommended that you use the default settings. Click "Next" when done.

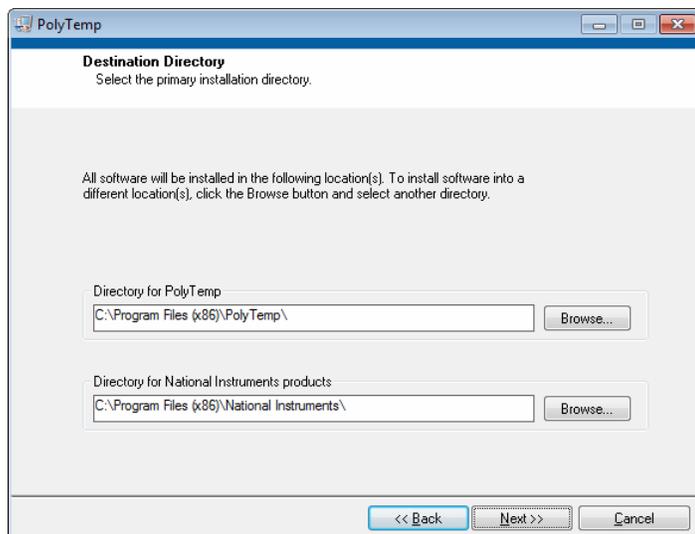

- 3) Select "I Accept the License Agreement" and click "Next" for both software licenses that appear. Note: If you choose not to accept the License Agreement, the installation will be cancelled.

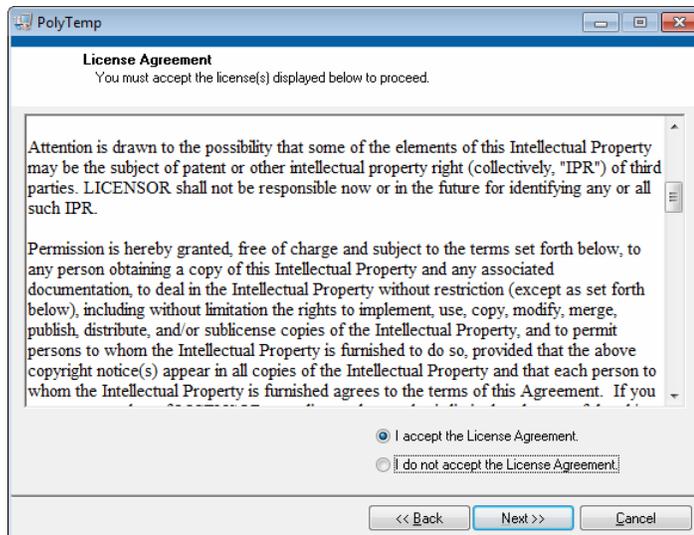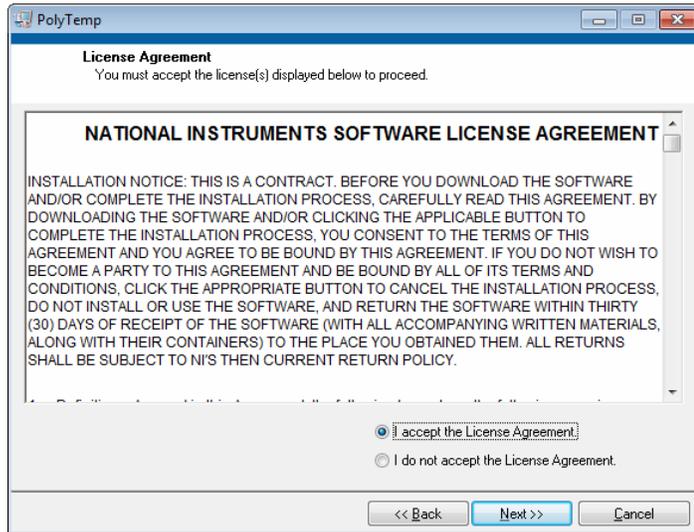


- 4) Click "Next" to start the installation. Click "Finish" when done. To finalize the installation, the PC must be restarted.

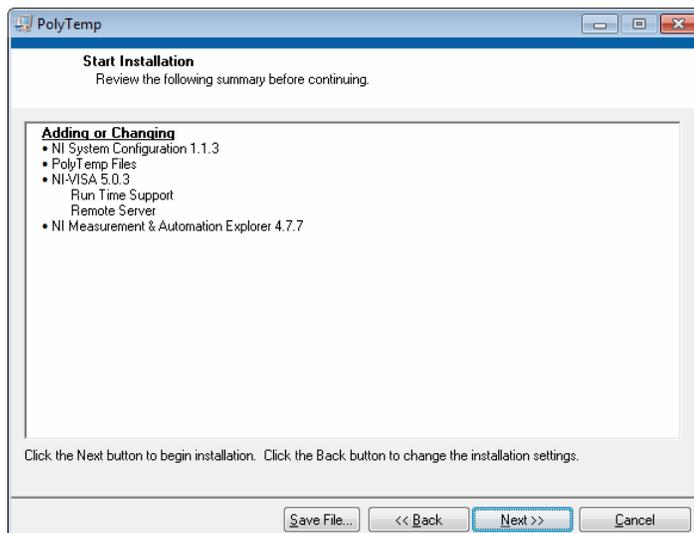

III. Installing the TDMS Viewer Plug-in for Microsoft Excel


IMPORTANT: PLEASE CLOSE ALL RUNNING APPLICATIONS ON YOUR PC BEFORE INSTALLING THE TDMS VIEWER.

- 1) Insert the PolyScience Driver CD into your PC.
- 2) Click the "Install TDMS Viewer" icon
- 3) Make sure all other applications are closed and click "Next"

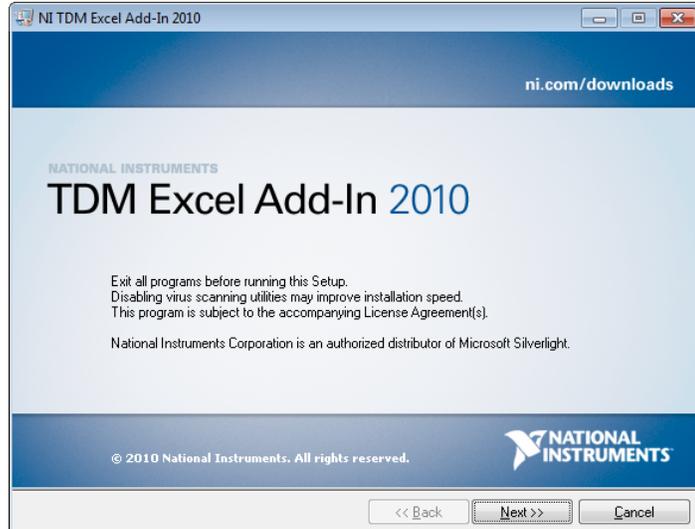


- 4) Select whether you would like to send data to National Instruments. NOTE: Unchecking this box will not affect the installation.

Click "Next" to proceed

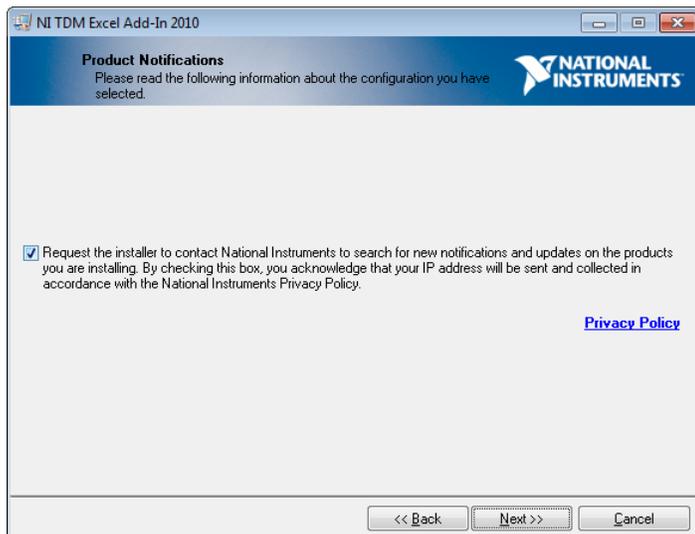

- 5) Select "I Accept the License Agreement" and click "Next" for both software licenses that appear. Note: If you choose not to accept the License Agreement, the installation will be cancelled.

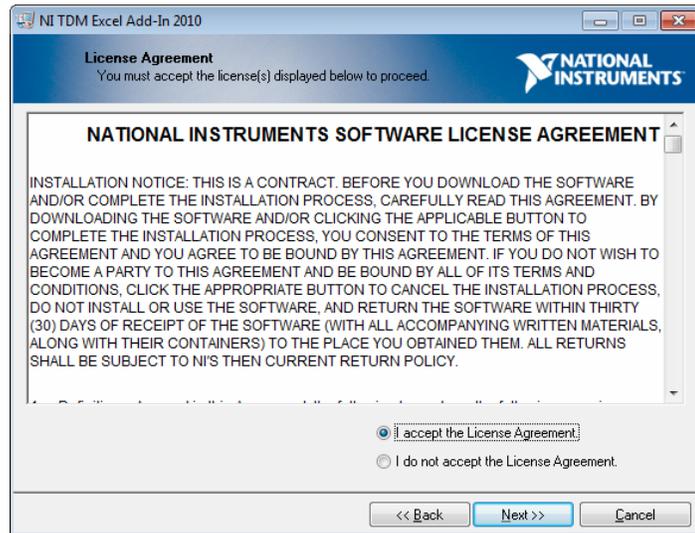

- 6) Click "Next" to start the installation. Click "Finish" when done.

